

PERFORM Operating Document

Use and Maintenance of QQQ LC-MS

PC-POD-CA-001-v04

Revision History

Version	Reason for Revision	Date
04	Condense version	06-Aug-2020

I. Introduction

The content of this PERFORM Operating Document (POD) provides guidelines for the use and maintenance of the QQQ LC-MS.

The PERFORM Centre Clinical Analysis suite has an Agilent 6460 Triple Quad Mass Spectrometer (LC-MS) which is a bench top QQQ providing atmospheric pressure ionization of liquid samples. A wide variety of compound classes may be analyzed using either ESI or APCI. Molecular weight information as well as structural information through collision induced dissociation provides useful qualitative sample information. This is well suited for both routine and complex analyses such as metabolism, toxicology, environmental, food safety, pharmaceutical, and drugs of abuse testing. This system provides analysis with greater sensitivity and productivity for quantifying pharmaceutical candidates, measuring trace-level environmental or food contaminants, or confirming biomarkers. Mass Hunter Quantitative Analysis software provides quantitative results with minimal manual intervention.

2. Definition of Terms and abbreviations

LC-MS	Liquid Chromatography-Mass Spectrometry
MS	Mass Spectrometer
QQQ	Triple Quadrupole
UPLC	Ultra Performance Liquid Chromatography
HPLC	High Performance Liquid Chromatography
N ₂	Nitrogen
ESI	Electrospray Ionization
APCI	Atmospheric Pressure Chemical Ionization
BinPump	Binary Pump
TCC	Thermostat Column Compartment
TFA	Trifluoroacetic acid
MP	Mobile Phase
EMV	Electron multiplier voltage
MW	Molecular weight
L	Liter
Min	Minute
mL	Milliliter
μA	Micro Ampere

3. Overview of LC-MS

- The I290 Infinity II LC System (UPLC I290) is equipped with a binary pump for two pairs of solvents (A1,A2 and B1, B2) and is capable to withstand 1200 Bars of operational pressure. Column and sample temperature is controlled by using column and auto-sampler thermostats. The chromatography system is coupled to mass spectrometer via ESI or APCI interfaces (sources) which are interchangeable.
- Two supplies of nitrogen are required to run MS analysis; One, from nitrogen tank (ultrapure grade) to operate the collision cell and another from N₂ generator to operate ESI or APCI source, The first source is equipped with filters which should be replaced at least once a year or more frequently if

required by workload. The second source of N₂ is supplied by a passive N₂ generator connected to a stationary source from Agilent. The pressure should be maintained sufficiently high to provide 80-100 psi pressure of nitrogen at the generator output. Diffusion/selection membrane in passive generator should be replaced at least once a year.

- There are 3 pumps (1 mechanical and 2 turbo) to create the vacuum in LC-MS system.
- There are 2 ion sources APCI and high efficiency Jet stream ion source ESI which are readily interchangeable.
- Built-in system logbook keeps log of errors for the sample running. Logbook deletes old entries each time a new worklist is run so it should be either saved or printed if required to report any malfunction of the instrument.
- Mass Hunter software suit is used for data analysis and is suitable for research and development (R&D) purposes. For compliance, additional software can be purchased.

3.1 Training requirements

Prior to using the LC-MS System individuals should:

- Read and sign this POD.
- Undergo appropriate LC-MS System training and/or provide a proof of an external training to the Supervisor, Clinical Analysis prior to use.

3.2 General precautions

- Use only water and high purity solvents (LC-MS grade): do not use water generated by Millipore system in the lab. Avoid contamination by using clean pipets, bottles and glassware.
- Make sure the level of oil in the mist filter from the mechanical pump is cleared regularly by opening the ballast valve for a few minutes. MS must be on standby before the ballast valve is opened. The valve should be closed for the routine operation. Make sure that there is no leak of oil. Check the oil level in the rough pump.
- If the rough pump overheats, it will beep. Let it cool by opening the front cover.
- The oil in the rough pump should be replaced and MS ion transfer capillary cleaned every 6 months. About 1.5 L of oil (Agilent 6040-0834) is required for replacement. To change the oil, turn off the LC-MS and shutdown the MS: In order and from the software, UPLC is on standby mode, right click on QQQ pane in the Data Acquisition software and press Vent. Wait 40 minutes. Then close the valve from the compress air (on the wall 90° clockwise). Turn off the computer, switch off the MS, UPLC, and the oil pump via their power buttons. The oil pump and capillary can be cleaned. To turn on the MS, open the compress air slowly and then manually turn on UPLC, MS and oil pump. Wait until UPLC injector finish the initialization. Turn on the computer, open Data acquisition software, right-click on Pump Down in the MS pane. Wait

PERFORM Centre

minimum 4 hours before using the MS and proceed to an autotune before any run.

- Verify the pressure from the Nitrogen generator, outlet pressure should not be below 80 psi, if the pressure is too low, the system will indicate an error and the system will not be able to operate the method.
- Check the calibrant level before tuning. Use different tuning mixes for both ESI and APCI sources available in the lab.
- Verify that all the cables are properly secured.
- It is important to use appropriate fittings with columns, valves and unions. Agilent Technologies uses Swagelok fittings but for Waters columns use Waters fittings.
- Capillary tubing is color coded to identify the internal diameter e.g. Red= 0.12 mm, Green= 0.17 mm and Blue = 0.25 mm. Make sure to use the right tubing and fittings for the connections.
- The binary pumps should be purged and primed before use.
- Drain any liquid from the LC-MS waste bottle otherwise it will flush back to LC-MS.
- Set the pressure limits according to your column.
Buffers should be <15 mM and sonicate prior to use. Never leave pure acetonitrile on the system as it will polymerize if left standing for periods of time. Always remove the column and store the system with 10% acetonitrile in HPLC grade water.
- MP bottles should never be covered with Para film. Use Aluminum foil.
- If the sample injection is required using a 96 well plates, configure the software for 96 well plates. By default it is configured for injection using vials.
- Flush the system with water right after if the buffer has been used to get rid of the salts. When changing solvents, all channels should be purged at 2 mL/min for 20 minutes/channel.
- If the column is not in use, rinse it with additive free mobile phase and then with 50-80 % acetonitrile at 0.1 mL /min for 10-20 column volumes or as recommended by the column manufacturer.
- If very sticky compounds are used, extra cleaning is required to remove traces. An Injector Purge Kit is available from Agilent Technologies for extra cleaning for sticky compounds.
- To change the APCI source put system on standby. Yellow light “should turn on” on standby mode. Wear gloves while replacing the regular spray shield for APCI because it may be hot.

3.3 Relevant documents

- Agilent 6460 Series LC-MS QQQ techniques and operation **volume 1** Student Manual
- Agilent 6460 Series LC-MS QQQ techniques and operation **volume 2** Student Manual
- Nitrogen generator Manual

4. Procedure

4.1 MS Operation

If the system is not turned on, follow the sequence below to turn it on:

- Check the pressure in N₂ generator first for the pressure above 80 psi.
- N₂ cylinder is always open for collision energy.
- Turn on the triple Quad.
- Turn on the UPLC and wait for 5 minutes.
- Login with the given user name and password.
- Go to the Mass Hunter Data Acquisition application and turn on the system.

Always perform auto tune once the system is turned “ON”.

If the mass spectrometer is not in use, keep in standby mode. Never turn the system off except for appropriate maintenance or cleaning.

Follow the instructions from page 195-204 of the student manual LC-MS QQQ techniques and operation volume I for preparing UPLC and sample preparation procedure. A QQQ routine pre-operation checklist is included in Appendix I and can be used to verify the LC-MS operation.

If any component of LC-MS system does not work inform the Supervisor, Clinical Analysis or delegate.

4.2 Data acquisition software layout and instrument controls

Open the data acquisition software. There is a default layout which can be used. The status of the instrument can be determined by looking at the instrument status pane. Green means that the instrument is ready for a run. Yellow means “not ready”, red means that an error has occurred in the module and you can check the logbook to view the error or hover over the red bar for an error description (inform the Clinical Analysis Suite supervisor). Blue means that the system is running and a pink color indicates that the instrument is waiting to inject.

Turn BinPump on and flush the injector. Ensure that lines are clear of bubbles, if not purge the lines via the BinPump pane. Flush the mobile phase from the binary pump for 60 sec. In TCC (thermostat column compartment), with direct injection, switch to column 1 or 2 depending on the choice of the column. Under “Bin Pump” icon you can set mobile phase composition, gradient steps and times, flow rate. Turn the pump on.

Save the data in R: drive with your folder. Contact PERFORM IT analyst in order to obtain access to this drive.

Refer to pages 365-418 of the student manual LC-MS QQQ techniques and

operation volume I to set up all UPLC and QQQ parameters.

Follow the procedure below if you need to perform a Checktune or Autotune on LC-MS.

4.3 Check Tune and Auto Tune

Refer to page 171 of student manual LC-MS QQQ techniques and operation volume I for Checktune and Autotune procedure.

- Wait at least 15 min to stabilize before tuning. There are different low concentrations tuning mix solutions for ESI and APCI sources from Agilent Technologies. The flow rate for tuning mix is 0.3 mL /min. Make sure that there is enough calibrant in the bottle for Checktune or Autotune.
- Check tune once a week when in use or as specified in the study protocol. If one or two masses do not pass then do Autotune.
- Autotune will automatically change the QQQ parameters and calibrate the QQQ. Routine use only requires resolution and mass axis verification for polarity and resolution modes to be used. Select the ions of interest as Positive or Negative for the Check tune. It will take 15-20 minutes for every mode. No need to inject anything as it is automatically run by the system. If there is no peak in the Autotune, a number of issues may be causing this: the source is dirty, ion transfer capillary is blocked, there's no calibrant solution left, and/or calibrant solution is not being delivered to the source. This is not an extensive list but can help to start troubleshooting. To clean source flush it with 50:50 methanol and HPLC grade water before repeating.
- After Autotune the EMV is usually 2000 V for old instrument. For the new instrument the EMV is around 1100 V.
- Check the results if they are acceptable or not. The report will be automatically saved as PDF under D:\Masshunter\tune\QQQ\G6460A.
- Autotune should be done monthly to optimize ion transmission, update EM voltage and minimize source exposure to TFA. It takes about 60 minutes to run the Autotune for both Positive and Negative modes. Flush source with LC flow after Autotune before running samples in negative ions.

If Checktune fails and abundances drop dramatically, in parallel with a major loss in assay sensitivity, this suggests a problem with the source and/or mass analyzer. A major drop in tuning mix mass without a loss in analyte signal suggests a calibrant problem. The calibrant level may be low or it has been degraded or the diverter valve may be malfunctioning. Inform Supervisor of Clinical Analysis Suite.

Note: It is useful to monitor the Autotune EMV in the Autotune report in order to know when the electron multiplier will need to be replaced. When the sum of the Autotune EMV and the method delta EMV approach 3000 volts, a new electron multiplier tube should be

replaced after venting the LC-MS system. Inform Supervisor, Clinical Analysis who will place a service call.

4.4 Work list

Refer to pages 5-27 of the training manual (volume 2) for the work list preparation.

4.5 Qualitative analysis

Refer to pages 52-104 of the training manual (volume 2) for the Qualitative data analysis preparation.

4.6 Quantitative Analysis

Refer to pages 178-240 of the training manual (volume 2) for the Qualitative data analysis preparation.

4.7 Parameters to use in the method optimization

For the method optimization, the parameters for Jet stream ESI and APCI can be found at: https://perform.concordia.ca/booking/Reference/clinical_analysis/start-up_parameters_for_qqq_LC-ms.pdf

4.8 Preventive maintenance

Follow **Agilent 6400 Maintenance Guide** for detailed procedures. In general, the following is required for the regular maintenance of LC-MS when in use:

Daily:

Flush the nebulizer, clean electrospray spray chamber, clean the APCI chamber, clean multimode source, check drain bottle.

Weekly:

Clean electrospray spray chamber, clean the APCI chamber, Clean multimode source, and check the level and color of the roughage pump fluid. Check the oil mist filter.

Monthly:

Flush the nebulizer. Replace the oil pump every six months. Replace it sooner if the oil appears dark or cloudy.

Annually:

Replace MS selection wall rotor seal. Multiplier needs replacement every year. Change Hydrocarbon filters. Wipe the skimmer once a year.

When required:

Clean and adjust electrospray needle as well as the APCI needle. Clean the

PERFORM Centre

capillary, clean skimmer, change lamp in the source, check for leaks, replace the LC filter elements, replace electron multiplier horn, replace high energy dynode. Add rough pump fluid when the pump fluid level is low. Replace the fuses. Change the seal of the pump when required, needle wash and needle seat after 60 000 injections.

To adjust the needle, put it in to the magnifier and adjust the needle so that it is aligned with the tip of ESI or APCI. If it is blocked, dip it in 50:50 methanol and Type I water and sonicate it for 3 minutes.

Nitrogen generator:

Refer to Nitrogen generator Manual for the operation, troubleshooting and maintenance details. Below is the maintenance schedule for N₂ generator:

Part	Action	Frequency
Filters	Replace carbon absorber	1x per year
Oxygen sensor	Replace oxygen sensor	1x per 3 years
Oxygen sensor	Calibrate oxygen sensor	1x per year
Air compressor	Replace on failure	Expected life > 8000 hrs
Nitrogen compressor	Replace on failure	Expected life > 8000 hrs
Battery on PCB	Replace on failure	Expected life > 5 years